


# Changing the silicone tube cut-off for Schilling burettes (Art.-no. 1 068, 1 069)

1. Use a knife to cut off the glass discharge tip.  
Remove any silicone remains carefully.
2. Remove the 3 screws at the Fitting with outlet tap.  
The burette part with the silicone tube can be removed.
3. Cut off the old tube at the glass olive of the burette, carefully remove the remains.
4. Push the replacement tube back onto the olive.
5. Open the stopcock system by pressing and push the glass part through. Silicone tube cut-off must not be squeezed.
6. Retighten the screws and reinsert the glass discharge tip into the silicone tube.

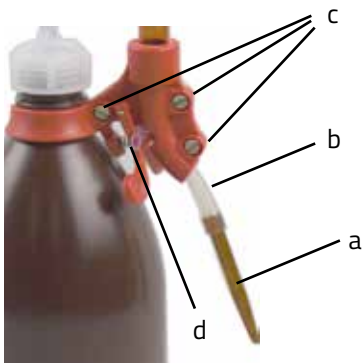

- a) Glass discharge tip
- b) silicone tube cut-off
- c) 3 screws
- d) PVC-tube



Spare Parts for Schilling Burettes		Art.-No.
Glass discharge tip for Schilling burettes (a)		1 075
Silicone tubes cut-off (b)		1 076
PVC-tubes cut-off (d)		1 078